

Update Instructions of the software qPCRsoftTOUCH V3.3.8.0 for qTOWER³/G touch

Download the setup data, unzip the folder or use the CD provided.

The download link is: <https://download.analytik-jena.de/?296a0e4c1a3c0b9f>

1. Switch on the qTOWER³touch and check the version of qPCRsoftTOUCH on the page *Options*. If the version is 3.3.8.0 no update is required.
2. Copy the folder *1-TabletPC* of the setup CD to a USB stick and insert it into the USB connector of the qTOWER³touch.
3. If the version is 3.3.4.0 quit qPCRsoftTOUCH by touching *Options* and then the exit button on the bottom right corner. The Windows desktop appears.
4. Find out which type of tablet PC is installed in your qTOWER³touch
 - a. **Device version 1 - (Odys /Trekstor):** With this type the tablet PC must be switched on separately on the right hand side of the upper hood.
 - b. **Device version 2 - (Avalue):** Available since 2019. With this type the tablet PC is switched on by the mains switch of the qTOWER³touch. On the right hand side of the hood only the USB connector is visible.

A: Update instruction for Device version 1 - (Odys / Trekstor)


- Use the Windows file explorer, select the USB drive, open the folder *1-TabletPC* and run the file *Setup_qPCRsoftTOUCH_V3-3-8-0.exe*. The following screen appears:



- Double touch the button *Device version 1 – (Odys/Trekstor)*.
- When you are asked for the administrator password enter “analytikjena”.
- Follow the instructions of the setup software.
- Note: You must run the setup twice because the first run uninstalls the old version only.
- Restart the tablet PC and qPCRsoftTOUCH will start with the updated version.

B: Update instruction for Device version 2 - (Avalue)

- Follow the instruction for *Device version 1*
- Hence, at the setup screen double touch *Device version 2 – (Avalue)*
- After the setup you must additionally run the app *SetKioskMode.exe* which displays the following screen:

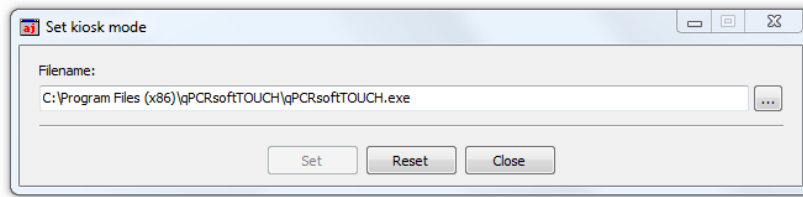

- Using the button [...] search for the file *qPCRsoftTOUCH.exe* in the Windows explorer. Usually it is found in the path shown in the figure above.
- Touch the button *Set* and note the message displayed.
- Restart the tablet PC and qPCRsoftTOUCH will be launched automatically in the updated version

The required app "*SetKioskMode.exe*" you can download here:

<https://download.analytik-jena.de/?9bdd683b189fbee1>

If you have questions, please contact service@analytik-jena.de